

Charles River Diagnostics through TAGCenter for non-Transnetyx users

NOTE: The captured screenshots on this page are representative of the steps below.

- 1** Place your order in the Charles River ordering platform, Laboratory Testing Management (LTM). Please follow instructions on criver.com/TAGCenter
- 2** Select "Sign In" on the TAGCenter then select "Skip".
- 3** Scan the barcode on the Charles River Cell Line Kit. Note: you have successfully scanned your Charles River Kit when you see the CRL**** barcode appear on the screen.
- 4** Drop your Charles River Cell Line Kit in the TAGCenter slot.
- 5** Confirm your Cell Line Kit was dropped via the confirmation screen.
- 6** Your Cell Line Kit will be shipped according to the TAGCenter schedule listed on the home screen. You will receive results via email.

